

ROWLEY
BIOCHEMICAL
INC.

DANVERS INDUSTRIAL PARK
10 ELECTRONICS AVENUE,
DANVERS MA 01923
TEL: 978-739-4883 FAX: 978-739-5640
www.rowleybio.com

L-761

EOSIN B

PROCEDURE:

1. Stain well-washed sections or slides, that have been stained with an alum hematoxylin, for 1-3 minutes in working Eosin B solution*. Dilute the working stain further if the stain is too intense.

* To prepare working stain, prepare just before use:

One part **Eosin B Stock Solution** L-761-1
Two parts distilled water .

If the diluted solution becomes cloudy, filter and add more stock solution..

2. Wash in several changes of water
3. Dehydrate, clear and mount.

Stain Results:

Nuclei	Blue
Cytoplasm	Pink
Collagen	Deep Pink
Smooth Muscle	Not quite as pink as collagen

References:

Clark, G.: Staining Procedures, Williams and Wilkins Company, Baltimore, 3rd Ed., C. 1973, p. 37.