ROWLEY BIOCHEMICAL INC.

DANVERS INDUSTRIAL PARK 10 ELECTRONICS AVENUE, DANVERS MA 01923

TEL: 978-739-4883 FAX: 978-739-5640

www.rowleybio.com

A-122 RAPID <u>GROCOTT'S METHENAMINE-SILVER NITRATE</u> <u>METHOD FOR FUNGI (GMS) AND PNUEMOCYSTIS CARINII</u> <u>(MODIFIED FOR MICROWAVE)</u>

FIXATION: Formalin, 10% Buffered Neutral
SECTION: Cut paraffin section at 4-6 microns

STAINING PROCEDURE:

- 1. Deparaffinize and hydrate to distilled water.
- 2. Oxidize in <u>Chromic Acid</u>, 10% (A-123-1), and microwave for 20 seconds at maximum power.
- 3. Let stand for 10 seconds in hot solution, then rinse in running tap water.
- 4. Place in Sodium Bisulfite, 1% Aq. (A-123-2), for 60 seconds.
- 5. Rinse in running tap water followed by 3 changes of distilled water.
- 6. Place slides in Methenamine Silver Nitrate Working Solution and microwave for 20 seconds at maximum power.

*To prepare working solution:

A. Methenamine – Silver Nitrate Stock Solution Silver Nitrate, 5%, Aq. (A-123-3) – 5ml Methenamine, 3%, Aq. (A-123-4) – 100ml

B. Methenamine Silver Nitrate Stock Working Solution

Methenamine-Silver Nitrate Stock Sol'n Stock 25 ml Distilled Water 25 ml **Sodium Borate 5%, Aq.** (A-123-5) 2 ml

Prepare fresh prior to use. Do not use if cloudv.

- 7. Let slides stand in hot solution which has been placed in a beaker of water at 60°C for 30 or more seconds. The slides should remain until they are pale brown.
- 8. Rinse slides in 3 changes of distilled water.
- 9. Tone in Gold Chloride 1%, aq. (SO-227) for 10 seconds.
- 10. Rinse slides in distilled water.
- 11. Place in Sodium Thiosulfate, 2% Aq. (A-121-6) for 1 minute.
- 12. Wash in running tap water for 2 minutes.
- 13. Place in Light Green Stock Solution, 0.2% (A-121-7), for 30 seconds.
- 14. Rinse in distilled water.
- 15. Dehydrate, clear and mount in resinous medium.

RESULTS:

Fungi	black, sharp detail
Pneumocystis carinii	black
Mucin	dark gray - taupe
Inner areas Mycelia, Hyphae	rose
Background	green, nale

REFERENCES:

Grocott, R.G.: Amer. J. Clin. Pathol., <u>25</u>:975 (1955).

AFIP Manual of Histologic Staining Methods; ed. G. Luna, 3rd.; New York:

McGraw-Hill Publications, c. 1968, p. 230

Reviewed and Approved

September 28 2023 • Supersedes All Other Dates