

ROWLEY
BIOCHEMICAL
INC.

DANVERS INDUSTRIAL PARK
10 ELECTRONICS AVENUE,
DANVERS MA 01923
TEL: 978-739-4883 FAX: 978-739-5640
www.rowleybio.com

L-759 **GILL'S HEMATOXYLIN #1**
FOR CYTOLOGY

FIXATION: Wet-fix in 95% Ethyl alcohol.

STAINING PROCEDURE:

1. Ethyl alcohol, 95% for 15 minutes.
2. Two changes distilled water, 10 dips each.
3. **Gill's Hematoxylin # 1** (L-759-1) for 2 minutes.
4. Two changes distilled water, 10 dips each.
5. **Scott's Tap Water Substitute** (L-759-4) for 1 minute.
6. Two changes distilled water, 10 dips each.
7. Ethyl alcohol, 95% 10 dips and counterstain.

RESULTS:

Chromatin Blue to Blue Black

REFERENCES:

Gill, G.W., Acta Cytol., 18:300-311, 1974